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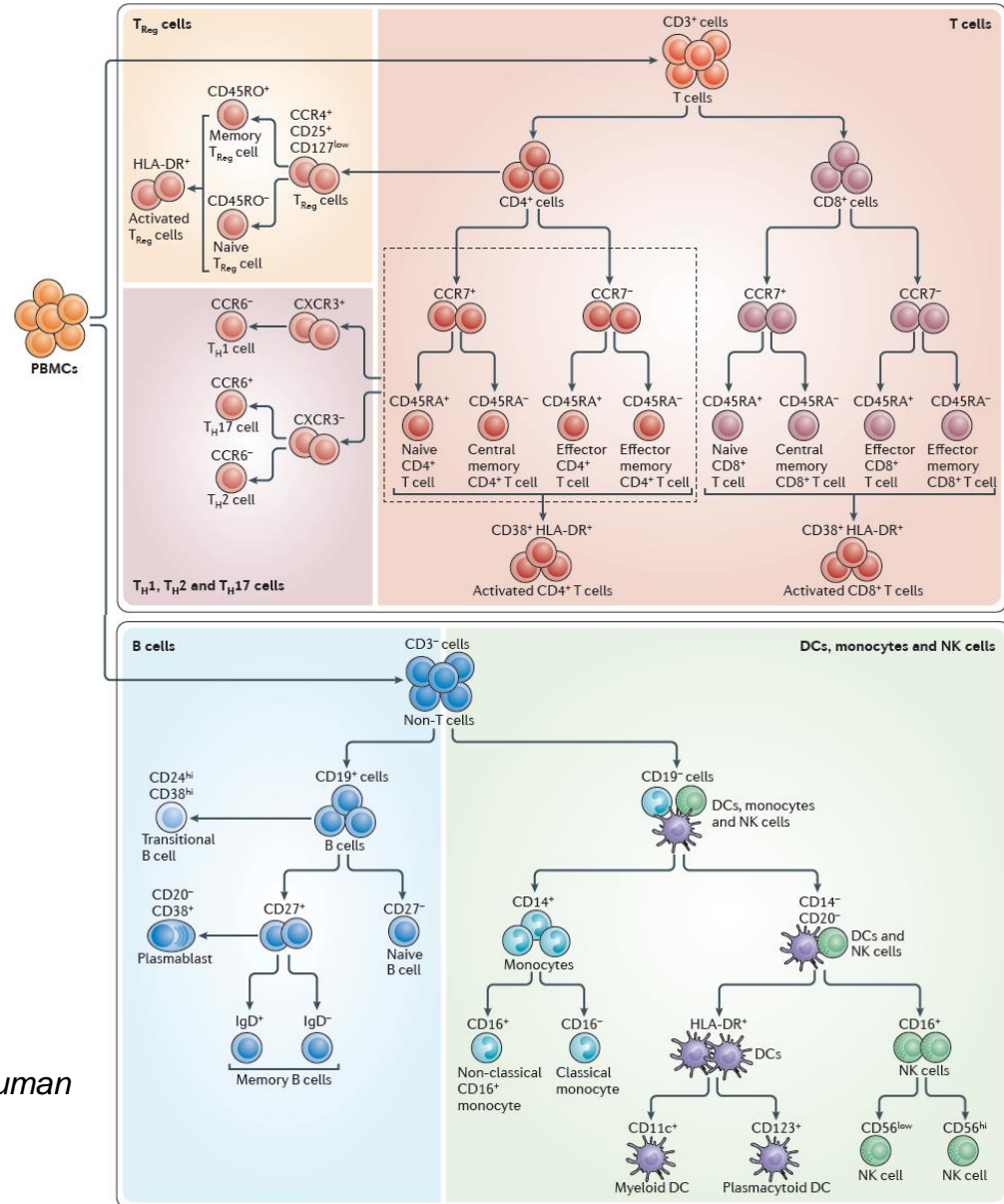
Standardizing immunophenotyping for the Human Immunology Project. Holden T et al.  
*Nature Rev Immunol* . ; 12(3): 191–200.

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
© 2016 Beckman Coulter, Inc. All rights reserved. Beckman Coulter, the stylized logo, and the Beckman Coulter product and service marks used herein are the trademarks or registered trademarks of Beckman Coulter, Inc. in the United States & other countries.

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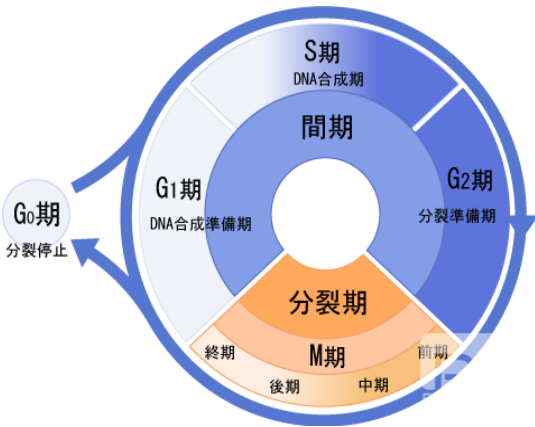
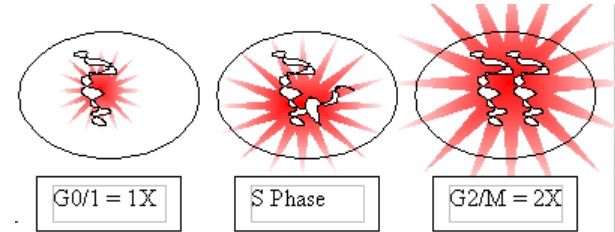
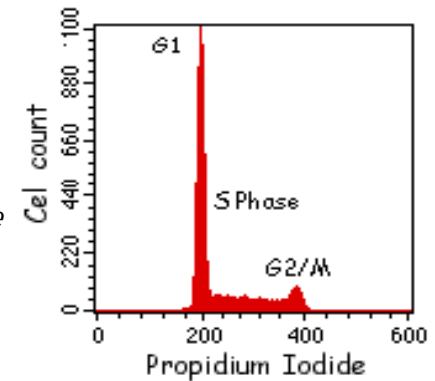
*These dyes are stoichiometric - number of bound molecules are equivalent to the number of DNA molecules present*

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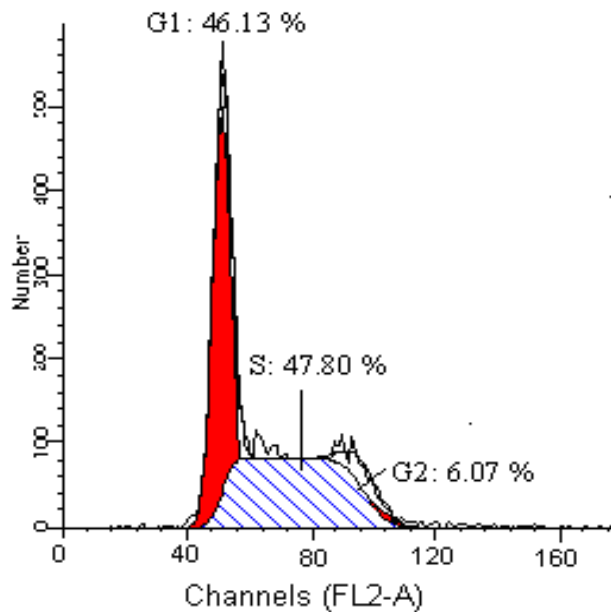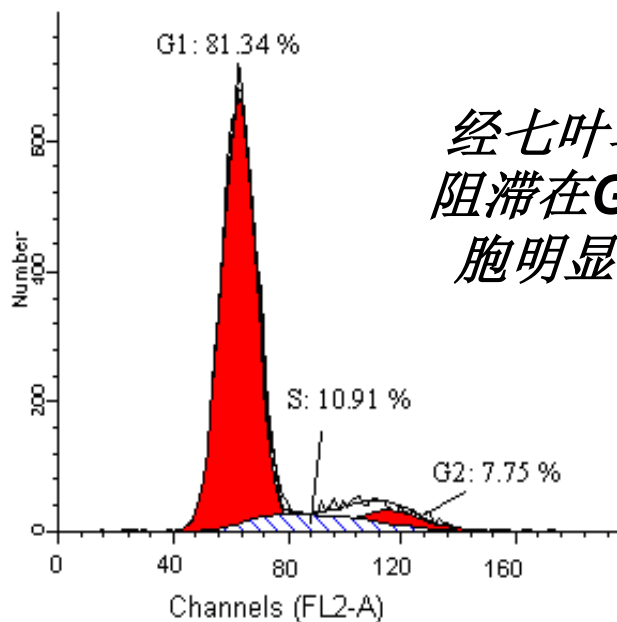
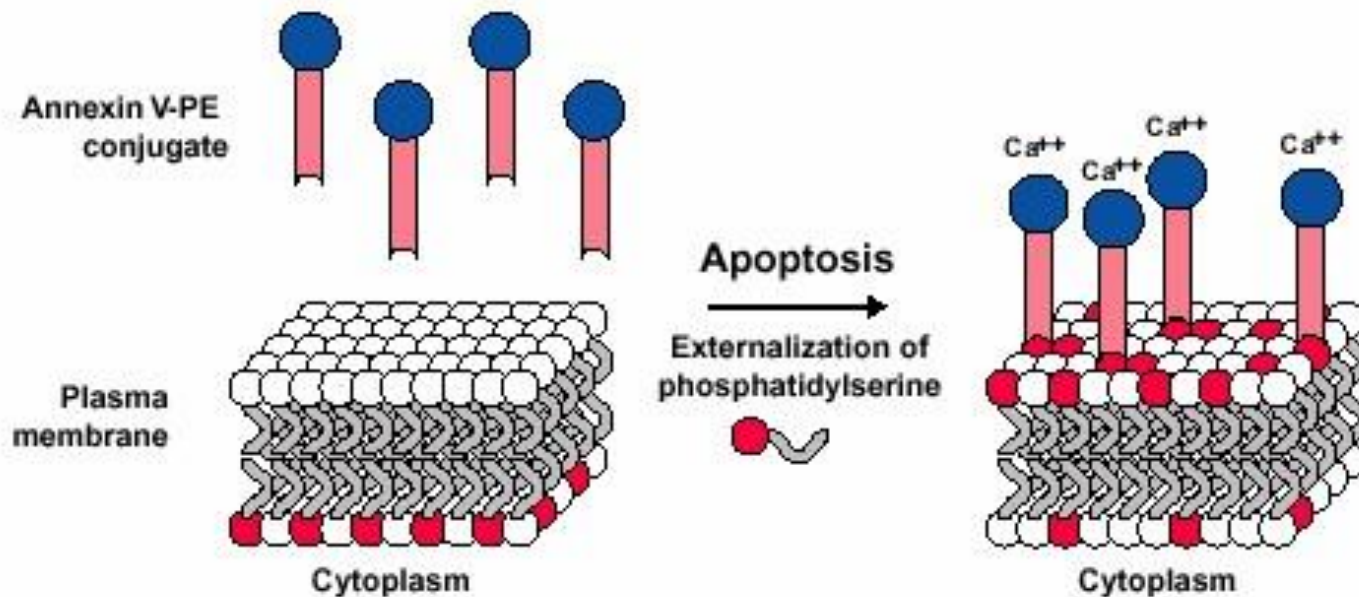
Schematic representation of the Annexin V assay.

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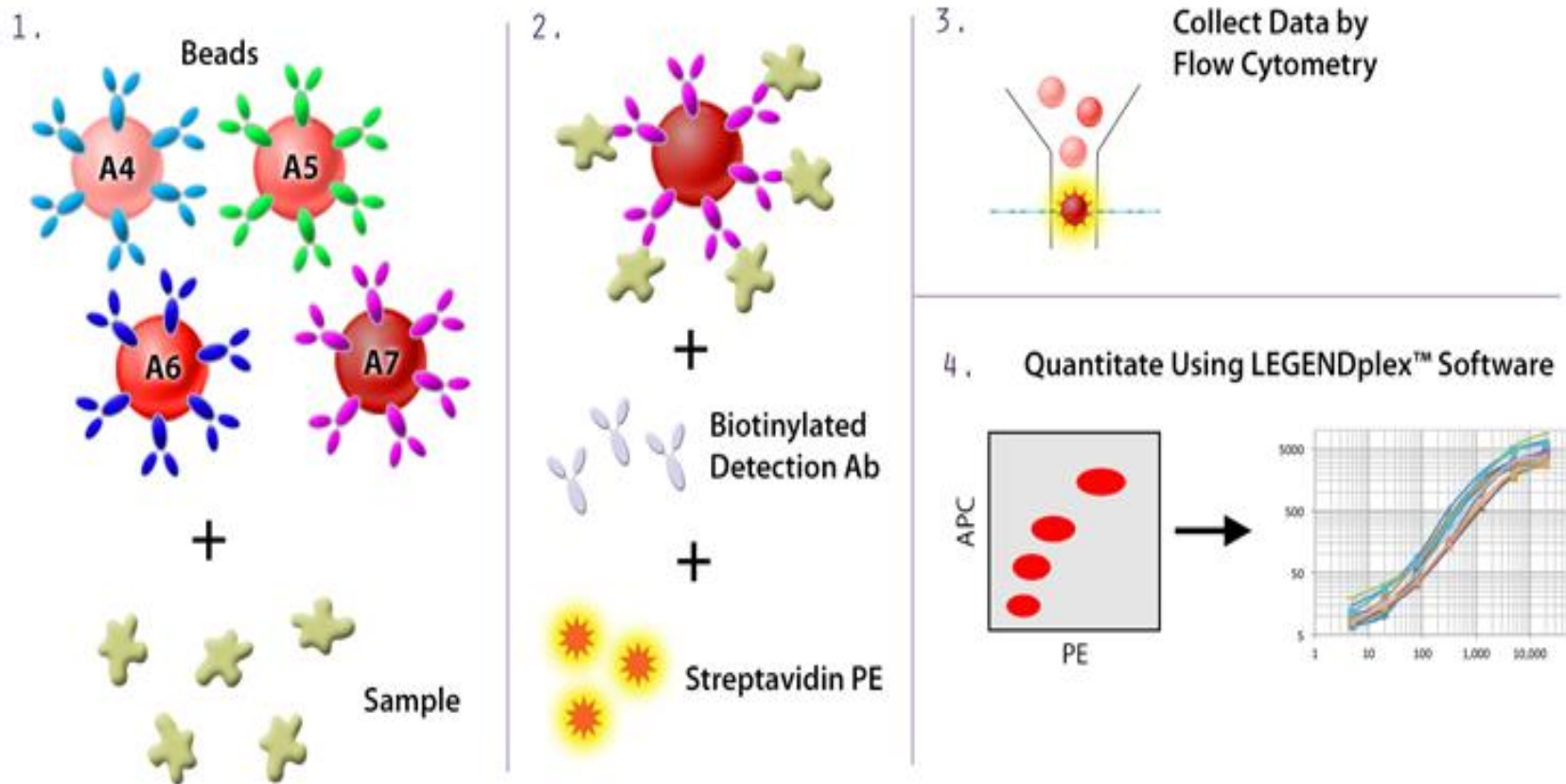
Delivering INNOVATIVE and trusted scientific solutions across the globe

©2016 Beckman Coulter, Inc. All rights reserved. Beckman Coulter, the stylized logo, and the Beckman Coulter product and service marks used herein are the trademarks or registered trademarks of Beckman Coulter, Inc. in the United States & other countries.

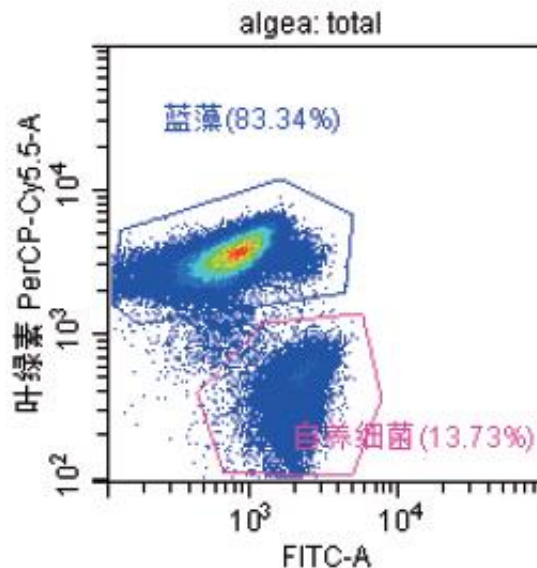
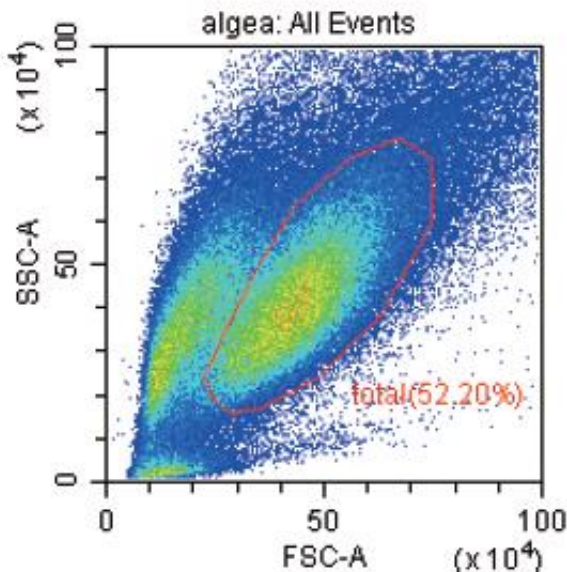
Internal / Company Confidential

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Tube Name: algea

Sample ID:

Volume(μL): 3.7

Population	Events	% Parent	Events/μL(V)
● 蓝藻	77090	83.34%	20704.57
● 自养细菌	12697	13.73%	3410.12

Delivering IT

©2016 Beckman Coulter, Inc. All rights reserved. Beckman Coulter, the stylized logo, and the Beckman Coulter product and service marks used herein are the trademarks or registered trademarks of Beckman Coulter, Inc. in the United States & other countries.

Confidential - Company Proprietary

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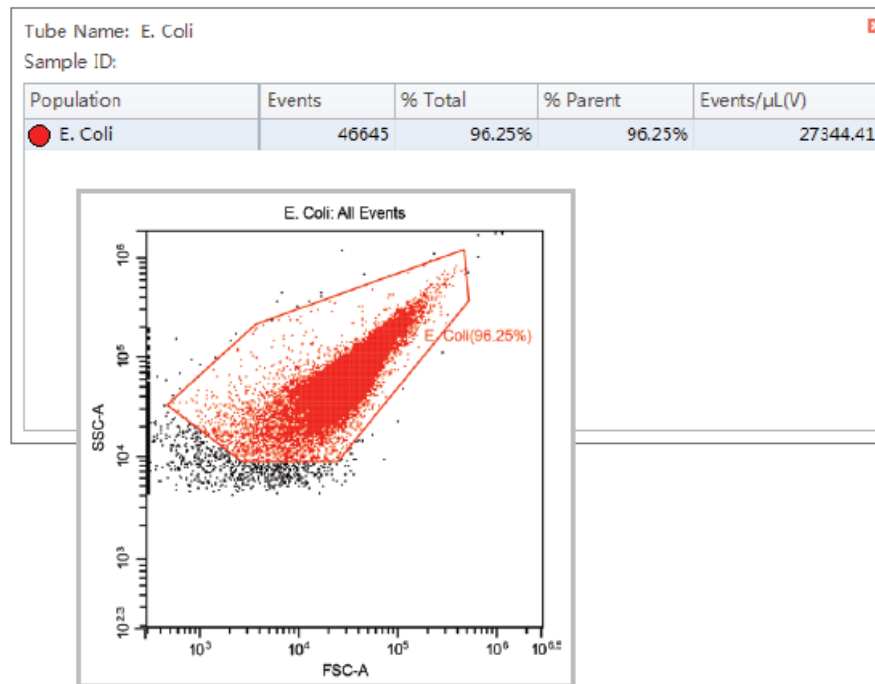
## Materials and tools

*E coli* sample

Delivering INNOVATIVE and trusted scientific solutions across the globe

©2016 Beckman Coulter, Inc. All rights reserved. Beckman Coulter, the stylized logo, and the Beckman Coulter product and service marks used herein are the trademarks or registered trademarks of Beckman Coulter, Inc. in the United States & other countries.

Confidential - Company Proprietary

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

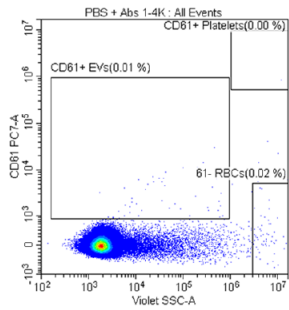
## A. PBS + Abs


## B. Whole Blood

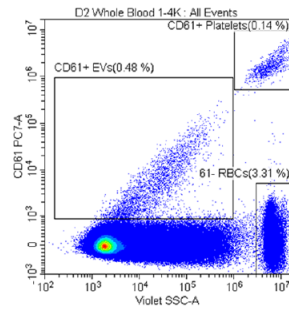
## C. Plasma

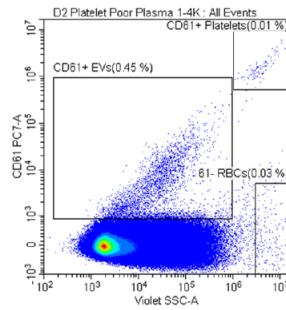
## D. Filtered Plasma

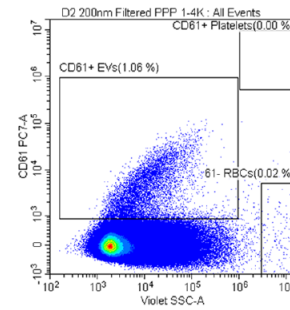
## E. Izon Fraction 5



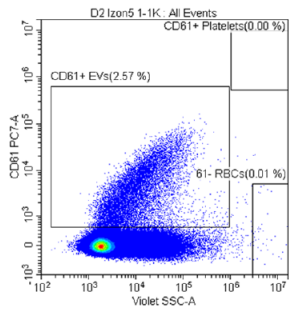
## F. Izon Fraction 6
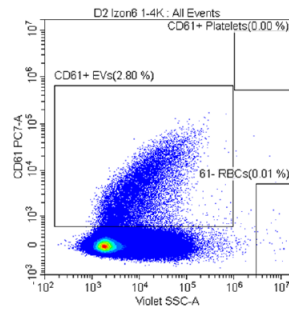
## G. Izon Fraction 7
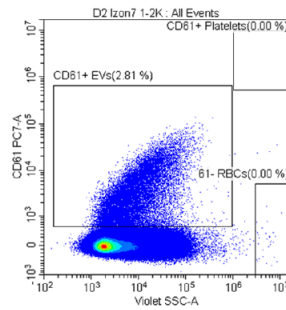
## H. Izon Fraction 8

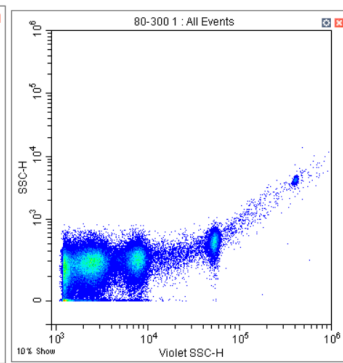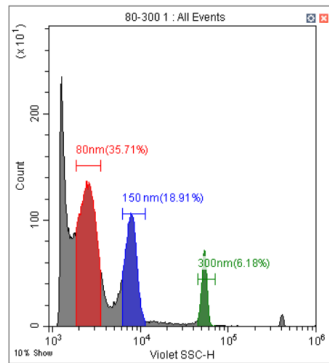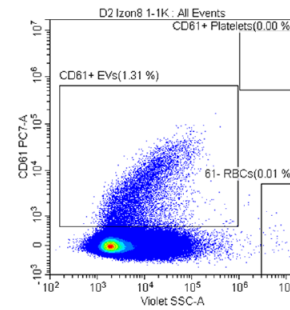Brittain, George C., et al. "A novel Semiconductor-Based flow cytometer with enhanced Light-Scatter Sensitivity for the Analysis of Biological nanoparticles." *Scientific reports* 9.1 (2019): 1-13.

Delivering INNOVATIVE and trusted scientific solutions across the globe

©2016 Beckman Coulter, Inc. All rights reserved. Beckman Coulter, the stylized logo, and the Beckman Coulter product and service marks used herein are the trademarks or registered trademarks of Beckman Coulter, Inc. in the United States & other countries.

Confidential - Company Proprietary

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

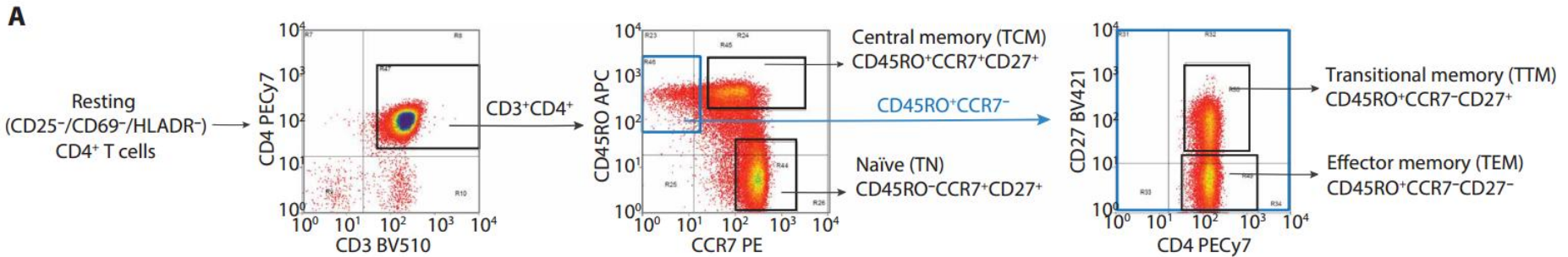
Different human resting memory CD4<sup>+</sup> T cell subsets show similar low inducibility of latent HIV-1 proviruses. Kwon KJ et al. Sci Transl Med. (2020)

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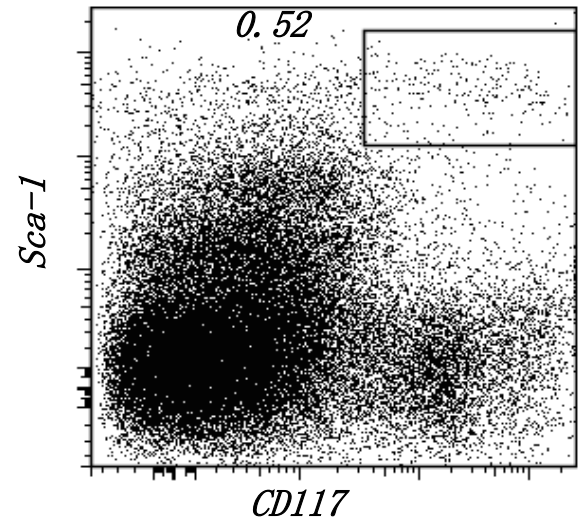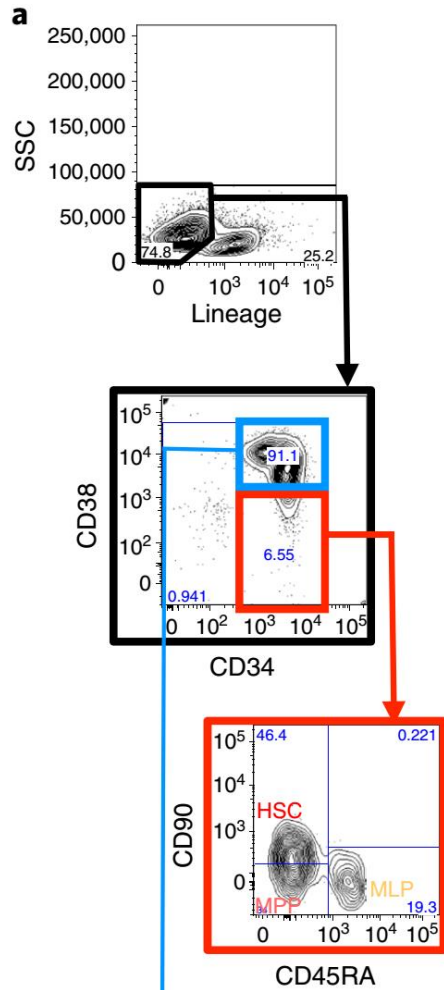
Dynamics of genetically engineered hematopoietic stem and progenitor cells after autologous transplantation in humans. Scala S et al. Nat Med. (2018)

Democratizing INNOVATIVE and trusted scientific solutions across the globe

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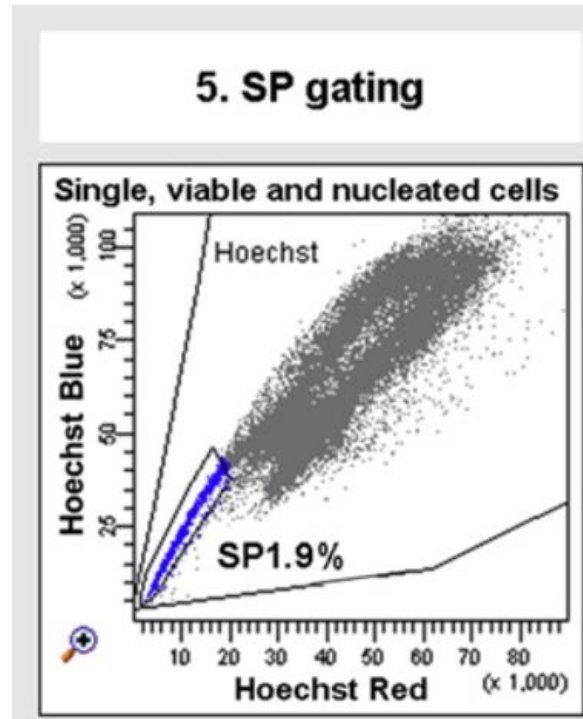Critical appraisal of the side population assay in stem cell and cancer stem cell research. Golebiewska A et al. Cell Stem Cell. (2011)

Delivering INNOVATIVE and trusted scientific solutions across the globe


©2016 Beckman Coulter, Inc. All rights reserved. Beckman Coulter, the stylized logo, and the Beckman Coulter product and service marks used herein are the trademarks or registered trademarks of Beckman Coulter, Inc. in the United States & other countries.

Confidential - Company Proprietary

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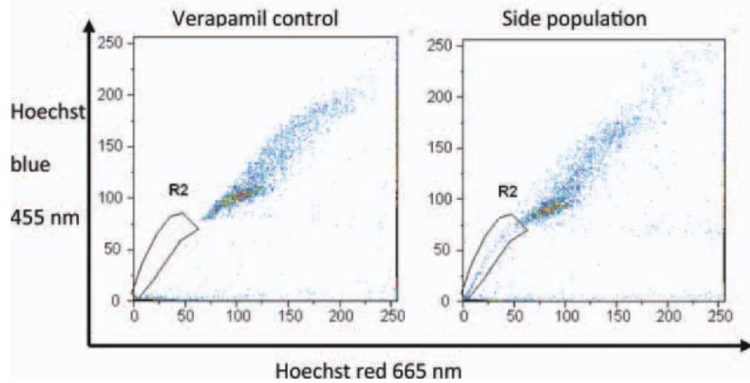**Table 1.** Overview of cancer stem cell markers suitable for flow cytometric analysis

	PUTATIVE MARKER PATTERNS FOR CANCER STEM CELLS (CSCS)	REFERENCES	ANNOTATION
Acute myelogenic leukemia (AML)	CD34 <sup>+</sup> /CD38 <sup>-</sup> CD90 <sup>+</sup>	(53,54) (55)	High risk AML
Acute lymphoblastic leukemia	CD34 <sup>+</sup> /CD38 <sup>-</sup> /CD19 <sup>+</sup>	(56)	
Bone sarcoma	Stro-1 <sup>+</sup> /CD105 <sup>+</sup> /CD44 <sup>+</sup>	(57)	
Brain tumor	CD133 <sup>+</sup>	(50,58)	
Breast cancer	ESA <sup>+</sup> /CD44 <sup>+</sup> /CD24 <sup>-low</sup> /lin <sup>-a</sup> )  CD90 <sup>low</sup> /CD44 <sup>+</sup>	(18)  (59)	Further enrichment by selection of ESA <sup>+</sup> cells Cells localized at the invasive front
Colon cancer	CD44 <sup>+</sup> /CD24 <sup>-low</sup> /ALDH1 <sup>high</sup> CD133 <sup>+</sup> ESA <sup>high</sup> /CD44 <sup>+</sup> /(CD166 <sup>+</sup> ) CD133 <sup>+</sup> /CD44 <sup>+</sup>	(51) (48,60) (61) (62,63)	
Colon cancer (metastatic)	CD133 <sup>+</sup> /CD24 <sup>+</sup> CD133 <sup>+</sup> /CD44 <sup>low</sup> /CD24 <sup>+</sup> CD133 <sup>-</sup> /CD44 <sup>+</sup> /CD24 <sup>-</sup>	(9) (64)	Two subsets of tumor-initiating cells identified
Endometrial cancer	CD133 <sup>+</sup> SP <sup>+</sup>	(65) (66)	
Gall bladder cancer	CD133 <sup>+</sup> /CD44 <sup>+</sup>	(67)	
Gastric cancer	CD44 <sup>+</sup>	(68)	
Liver cancer	CD133 <sup>+</sup> /CD44 <sup>+</sup> CD90 <sup>+</sup> EpCAM <sup>+</sup> CD133 <sup>+</sup>	(69) (70) (71) (24,72)	
Metastatic melanoma	CD20 <sup>+</sup>	(34)	
Ovarian cancer	CD133 <sup>+</sup> /ALDH <sup>+</sup> CD44 <sup>+</sup> /CD117 <sup>+</sup>	(73) (74)	
Pancreatic cancer	CD44 <sup>+</sup> /CD24 <sup>+</sup> /ESA <sup>+</sup>	(75)	
Prostate cancer	CD44 <sup>+</sup> /α <sub>2</sub> β <sub>1</sub> <sup>hi</sup> /CD133 <sup>+</sup> CD44 <sup>+</sup> /CD24 <sup>-</sup> SP <sup>+</sup>	(20) (76) (77)	
Renal cell carcinoma	CD105 <sup>+</sup> /(CD133 <sup>-</sup> /CD24 <sup>-</sup> )	(78)	
Head and neck cancer	CD44 <sup>+</sup>	(79)	No further enrichment by endothelial-specific antigen or CD24




<sup>a</sup> lin = lineage marker CD3, CD14, CD16, CD19, CD20, and CD56.

Flow cytometry in cancer stem cell analysis and separation. Greve B et al. Cytometry A. (2012)

Delivering INNOVATIVE and trusted scientific solutions across the globe

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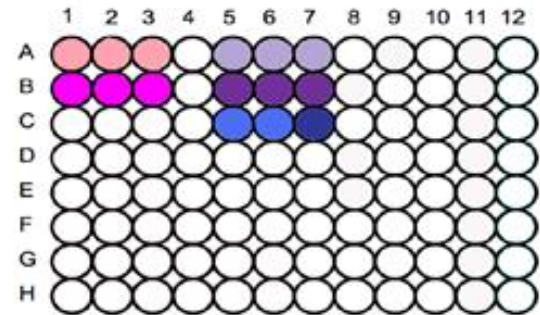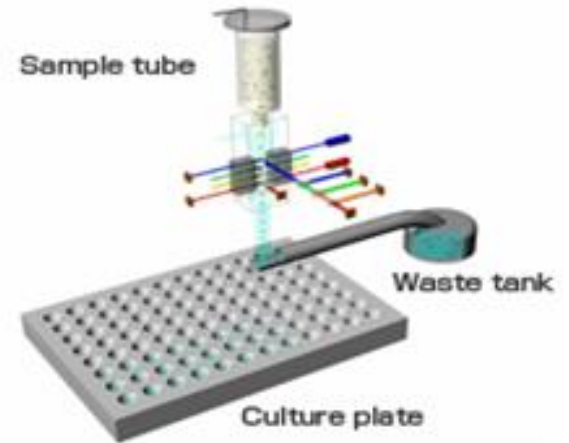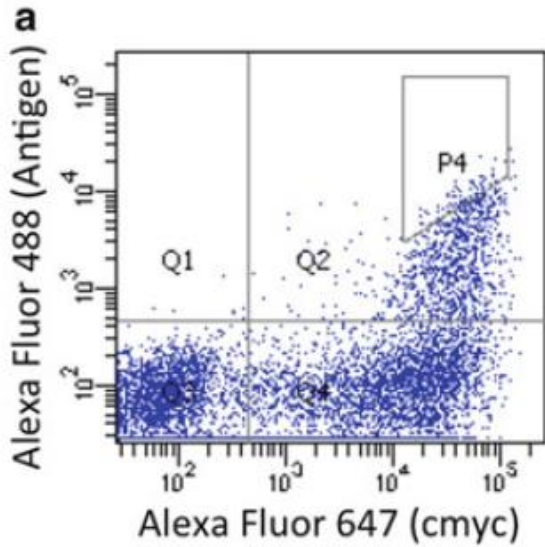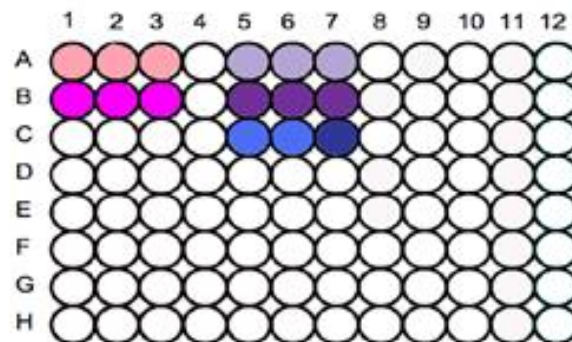
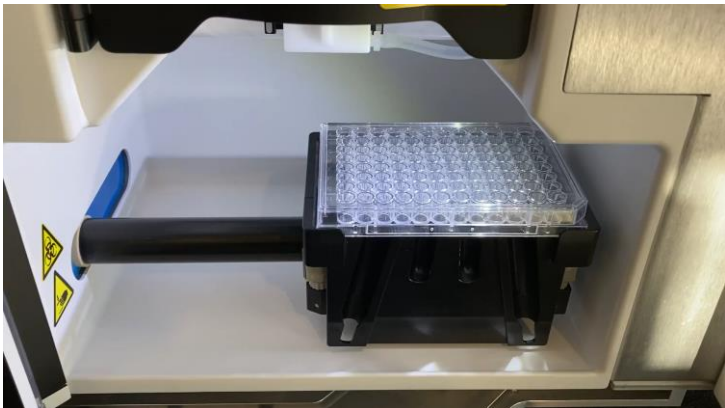
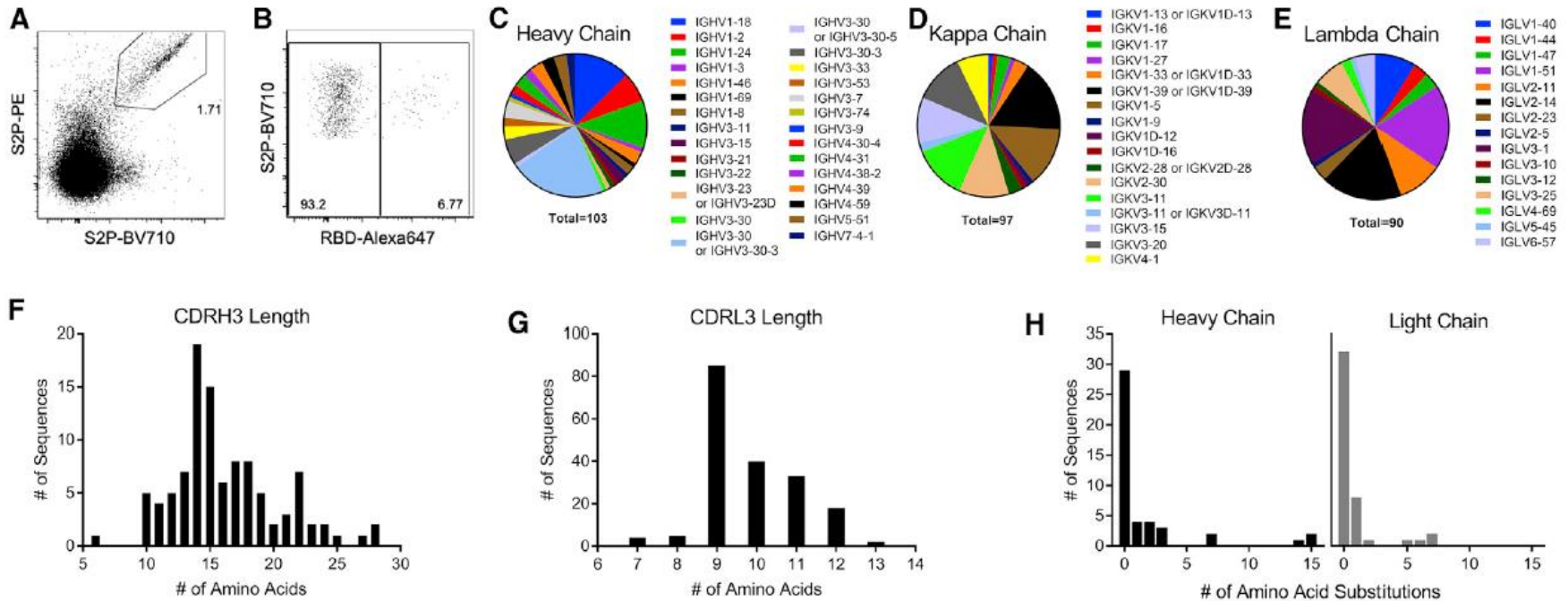
Figure 2. Early B Cell Response to SARS-CoV-2 Is Diverse and Largely Unmutated

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RESEARCH

Open Access

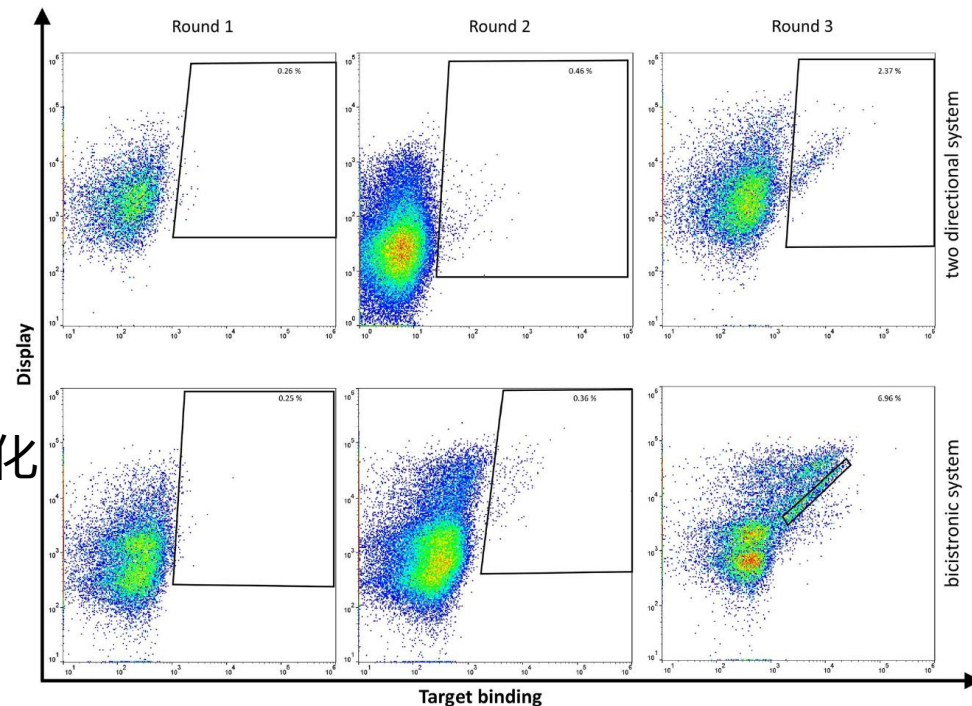
## A novel one-step approach for the construction of yeast surface display Fab antibody libraries

Simon Rosowski<sup>1</sup>, Stefan Becker<sup>2</sup>, Lars Toleikis<sup>2</sup>, Bernhard Valldorf<sup>2</sup>, Julius Grzeschik<sup>1</sup>, Deniz Demir<sup>2</sup>, Iris Willenbücher<sup>2</sup>, Ramona Gaa<sup>2</sup>, Harald Kolmar<sup>1</sup>, Stefan Zielonka<sup>2\*</sup> and Simon Krahl<sup>2\*</sup>

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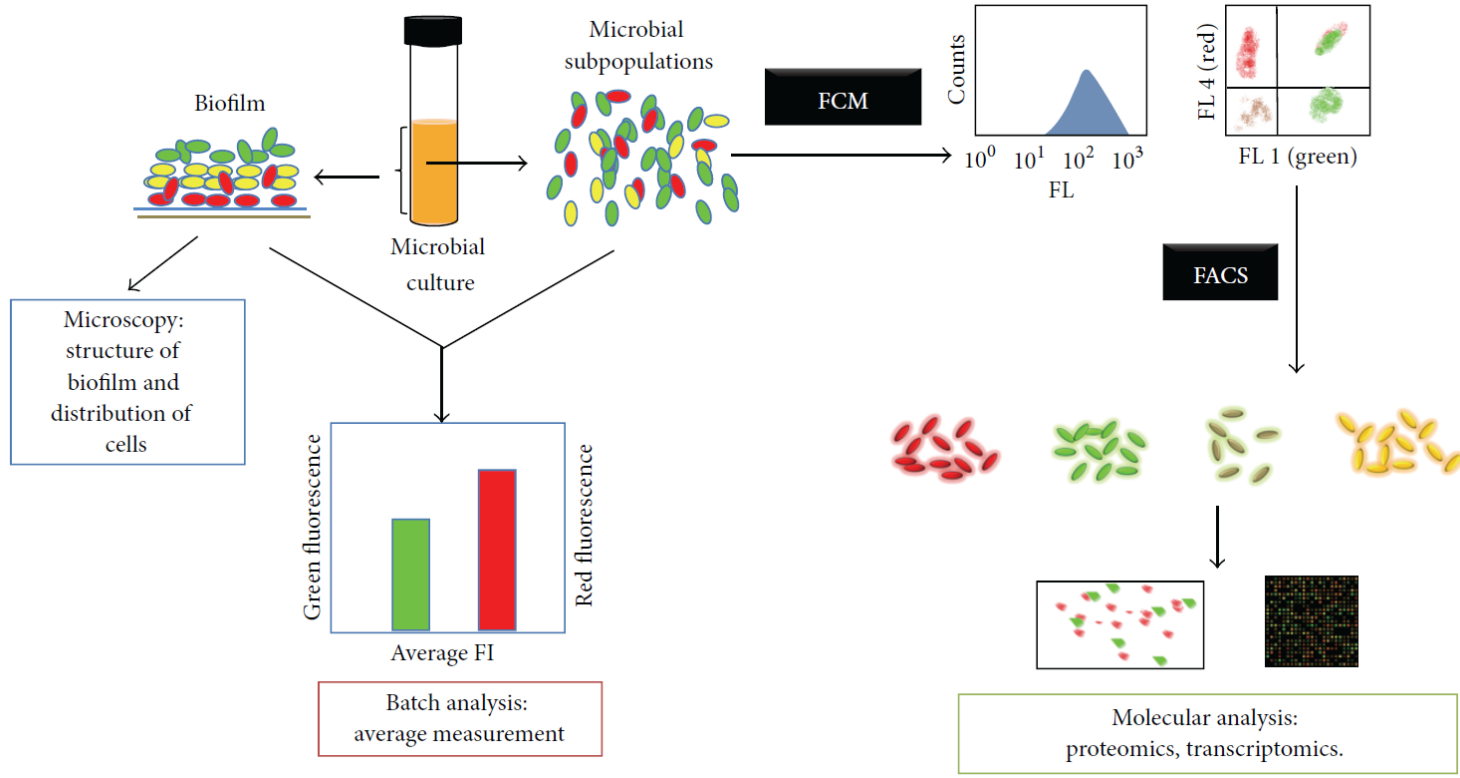
FCM allows analyzing subpopulations of cultured bacterial cells with different physiological states

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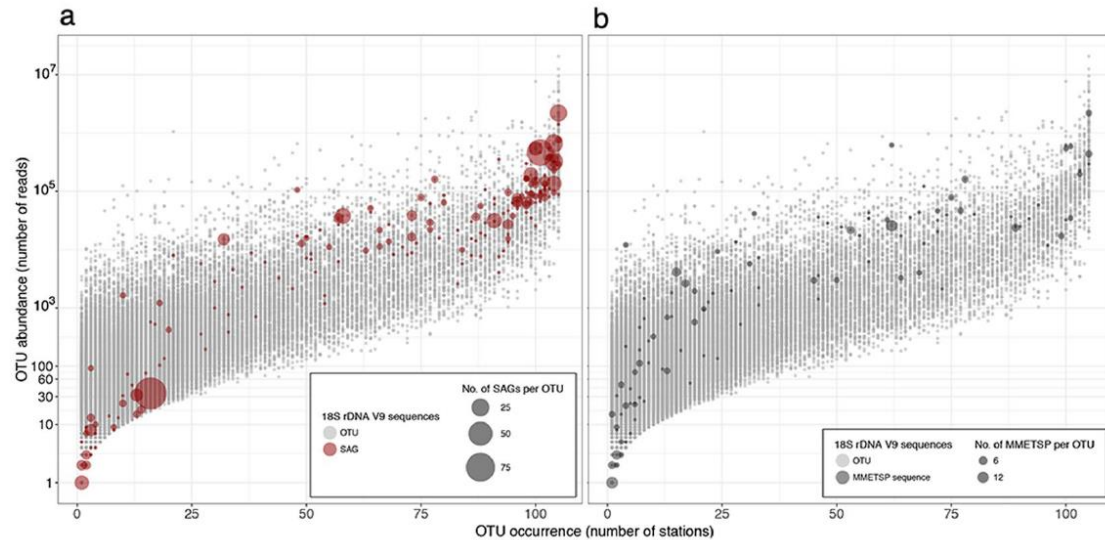
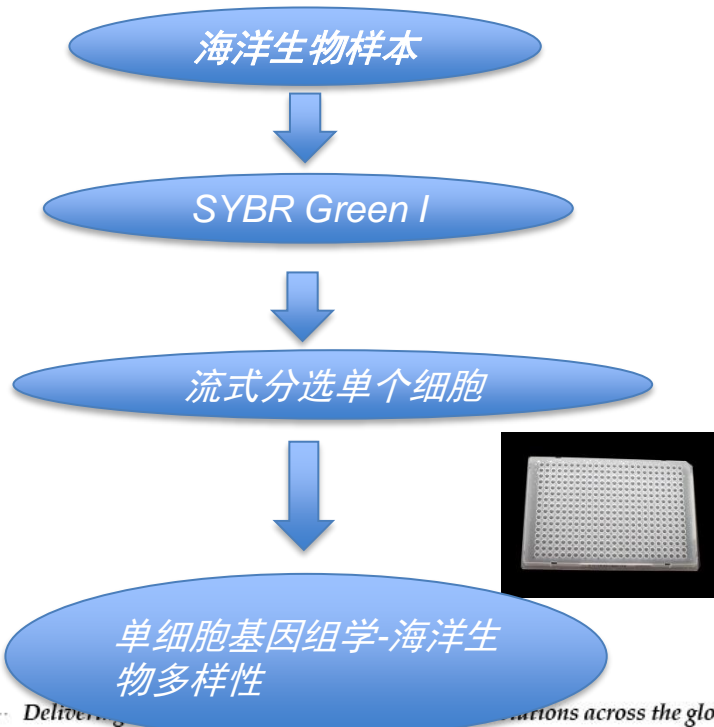
SCIENTIFIC REPORTS

OPEN Single cell genomics yields a wide diversity of small planktonic protists across major ocean ecosystems

vol. 1 June 2018  
doi:10.1038/s41598-018-28303-9

Delivering... stations across the globe

© 2016 Beckman Coulter, Inc. All Rights Reserved. Beckman Coulter, the stylized logo, and the Beckman Coulter product and service marks used herein are the trademarks or registered trademarks of Beckman Coulter, Inc. in the United States & other countries.

Confidential - Company Proprietary

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